

Suggested PCR Protocol

Pipette the following into a PCR tube, mix and make up to a final volume of 50 µL. We recommend dispensing all reagents on ice, adding the enzyme last. It is important to vortex all buffers and MgCl₂ solutions before use to remove any gradients that may result from repeated freeze/thaw steps.

If you do have already your own PCR-Protocol established, please use your existing pipetting scheme and Thermocycler protocol.

Components	Quantities
Template DNA	1 ng - 10 ng plasmid DNA or 5 ng - 500 ng genomic DNA
Nucleotides	1 µL (10 mM) each dNTP
10X amplification buffer	5 µL
primer 1:	4-7 µL of 3 µM solution (10 - 20 pmole absolute)
primer 2:	4-7 µL of 3 µM solution (10 - 20 pmole absolute)
sterile, bidistilled water	up to 50 µL
<i>ReproFast</i> -Polymerase	0.25 - 0.5 µL (1.25 - 2.5 units)

The following cycling protocol has to be seen as suggestion. Cycle numbers and cycle times have to be adjusted.

Cycling Profile *	
94°C for 5 minutes	step 1
94°C for 30 sec.	1 - 30 cycle(s)
annealing temp. for 15 sec.	1 - 30 cycle(s)
72°C for 15-30 sec. **	1 - 30 cycle(s)
94°C for 30 sec.	
72°C for 5-10 minutes	last step
cool down to 4°C	

* cycling times are proposals and have to be adjusted to the special needs.

** the longer the template to be amplified, the longer the amplification time

After running the amplification reaction, 5 µL of each sample should be analyzed in a 1% agarose gel containing 0.5 µg/mL of ethidium bromide.

Note: For every template/primer pair the optimal reaction conditions have to be evaluated empirically, changing the primer/template ratio, the ionic strength (with MgSO₄) and the cycle parameters (time and temperatures).

Trouble shooting:

No PCR® Product

Check if annealing temperature has been calculated correctly. The optimal annealing temperature of primers lies about 5°C below the T_m of the oligo.

TM calculation: For oligos up to 15 bases: $TM = 4x(G+C) + 2x(A+T)$
 For larger oligos: $TM = 81.5 + 16.6([\log_{10}J^*]) + 0.4(\%G+C) - (600/n) - 0.63(\%FA)$

A = Adenosine, C = Cytosine, G = Guanine, T = Thymidine, FA = Formamide, J* = concentration of monovalent cations, n = number of bases; T_m = calculated melting temperature,

Remark: If primers with mismatches are used the *ReproFast* Polymerase **must** be added immediately before the start of the PCR®. Do **not** let the reaction mix stand at room temperature since primers will be degraded by the *ReproFast* Polymerase.

Bands smear over the Gel

Increase annealing temperature

Add up to 5% DMSO to the Reaction mixture

Reduce the amount of template DNA

Check if the primers bind more than once on the template DNA. Perform control reactions with only one of the primers.

Low Yields

Increase annealing temperature

The PCR process is licensed and covered under US-Patent owned by Hoffmann-La Roche, Roche Int. USA